



Flexible and versatile design for optimized staining results

Devoted to Histology



## AUTOMATED SLIDE STAINER

# **SS-30**

### FOR ROUTINE AND SPECIAL STAINING PROTOCOLS IN HISTOLOGY AND CYTOLOGY LABS

Flexible, optimized and user-friendly

#### MULTISTAINING CAPABILITY

More staining flexibility allowing simultaneous and automatic staining of various 30-slide racks with identical or different staining protocols. Continuous loading to maximize efficiency and productivity in labs.

Up to 5 slide racks can be run simultaneously depending on protocols, load frequency and instrument configuration, enabling up to 150 slides to be stained simultaneously. The instrument calculates the most efficient route for each rack/ protocol.

#### FLEXIBLE STAINING

The agitation system can be individually programmed for each station: it is available in 4 different preset dip modes with 3 configurable parameters: number, speed and amplitude of the dips. A special programming for the washing stations is also available.

Several adaptors are available for the most popular coverslipper racks in the market.

#### COMPACT FOOTPRINT

Ideal for labs with shallow benches to optimize valuable space. Smartly designed to occupy less space without compromising slide staining productivity.

#### **REAL-TIME DISPLAY**

SS-30 displays in real time the status of each staining protocol. It can store up to 20 programmed protocols in memory with 50 steps each maximum as well as up to 52 reagents (32 pre-programmed).

The drain function minimizes the reagent carryover and helps reducing reagent consumption to optimize costs.

#### DEDICATED SOFTWARE

SS-30 stainer is supplied with a PC dedicated software that offers the option to define protocols and modify the list of saved reagents from the convenience of a PC.

It has also reporting capabilities through a USB plug to acquire data, displaying real status of the running protocols and managing reports and errors.



# **USER AND SPECIMEN SAFETY**

Lab-friendly integrated filtration system equipped with a charcoal filter and individual lids for the reagent stations to avoid evaporation of reagents. Optional adaptor for the laboratory general extraction system.

Back-up battery with 2h autonomy to finish the staining protocols in case of power failure. External 12V power supply.





The transparent hood is equipped with two hinged doors (loading/ unloading stations) to reduce the exposure to reagents.

Water overflowing in the stainer is secured thanks to an innovative two water level sensors system.

The Reagent Management System (RMS) keeps track of reagent usage and ensures an optimal staining quality (available only in version SS-30H).







#### **SPECIFICATIONS**

Specimens	Histological and/or cytological		
Staining capacity	Up t	o 5 racks simu	ultaneously
Slide rack capacity			30 slides
Number of programs		Up to 20 pro max. 50	grams with steps each
Immersion time	From	1 s to 59 m 59	) s per step
Agitation system	Indep	endently prog for e	grammable ach station
Selectable parameters			speed and ude of dips
Drain function	Min	imizes reagen	t carryover
Total stations	20 \	vessels with in	dividual lid
Reagent stations			Up to 19
Reagent stations capacity	/		300 ml
Water stations			Up to 3
Load stations			Up to 2
Unloading stations	(2 if d	rying station i	Up to 3 s available)
Drying station	(	available only	1 in SS-30H)
Drying station temperatu	ire		30 to 70° C
Fume extraction system		Cha	arcoal filter
Power requirements SS-30 SS-30H	100-240VAC 100-240VAC	50-60Hz 50-60Hz	0,17A 0,76A
Backup battery capacity			2 h
Dimensions (l x w x h)	1.200 x 440 x 368 mm		
Weight			55 kg

#### **ORDERING INFORMATION**

SS-30	Automated slide stainer	100-240v / 50-60 Hz
SS-30H	Automated slide stainer	100-240v / 50-60 Hz
	with drying station	

#### **OPTIONAL ACCESSORIES**

SS30-059	Adapter for THERMO-MICROM rack (30 slides)
SS30-200	Load station LEICA racks
SS30-200B	Lid for load station LEICA rack
SS30-201	Adapter for LEICA plastic rack ( 30 slides)
SS30-201B	Adapter for LEICA metal rack (30 slides)
SS30-202	Adapter for SAKURA rack (20 slides)
SS30-204	Adapter for MEDITE rack (20 slides)
SS30-205	Adapter for MEDITE rack (30 slides)
SS30-210	Megaslides adapter

#### Especialidades Médicas Myr S.L

ISO 9001 / 13485 certified company Lleida, 17-23 43712 Llorenç del Penedès - Spain

Tel. +34 977 66 8020 Fax. +34 977 66 8030 esp.medicas@myr.com.es www.myr.com.es

Local Distributor. Contact information.